

Corning® Hepatocyte Maintenance Medium for Cell Plating, Induction, and Hepatotoxicity Studies

Assay Protocol

CORNING



Materials

- ▶ Corning hepatocyte maintenance medium (Corning Cat. No. 40-550-CV)
- ▶ Cryogenic preserved primary human hepatocytes
- ▶ Cryogenic Preserved Hepatocyte Recovery Medium (CHRM)
- ▶ Collagen I-coated plates (Plate format options: 6-well plate, Corning Cat. No. 356400; 12-well plate, Corning Cat. No. 354500; 24-well plate, Corning Cat. No. 354408, and 96-well microplate, Corning Cat. No. 354407)
- ▶ Corning® Matrigel® matrix (Corning Cat. No. 356237)
- ▶ Multi-channel pipettors and tips
- ▶ Vi-CELL® cell viability analyzer (Beckman Coulter)
- ▶ Cell culture incubator (5% CO₂, 37°C, 95% humidity)
- ▶ Certified biosafety cabinet

Procedure for plating primary human hepatocytes

Day 1 (cell plating)

- Step 1:** Prepare hepatocyte plating medium by supplementing Corning hepatocyte maintenance medium with 10% FBS.
- Step 2:** Thaw a vial of cryogenic preserved primary human hepatocytes using CHRM medium according to the manufacturer's protocol.
- Step 3:** Re-suspend the resulting hepatocyte pellet in plating medium prepared in step 1.
- Step 4:** Estimate the density and percentage of viable cells by the trypan blue exclusion method using the Vi-CELL analyzer.
- Step 5:** Dilute the cells from previous steps with plating medium to a desired seeding density.
- Step 6:** Pipet cells onto a Corning Collagen I-coated plate(s) according to Table 1.

Plate	Viable Cells/Well	Volume/Well
6-well	1.4 x 10 ⁶	2.0 mL
12-well	0.7 x 10 ⁶	1.0 mL
24-well	0.35 x 10 ⁶	0.5 mL
96-well	0.06 x 10 ⁶	0.1 mL

Table 1. Recommended densities for plating hepatocytes

- Step 7:** Transfer the plates to the cell culture incubator and allow cells to attach to the plates for 6 hours.
- Step 8:** At the end of 6 hours, replace the plating medium with pre-warmed Corning hepatocyte maintenance medium.
- Step 9:** Optional Matrigel matrix overlay method: at the end of 6 hours, add the same volume of cold medium (approximately 4°C) containing 0.25 mg/mL Matrigel matrix to each well.
- Step 10:** Return the plates to the incubator for further experiments as required.

Procedure for CYP induction studies

Day 2 to 4 (3-day induction)

Step 11: Remove the culture medium from each well and replace with pre-warmed Corning® hepatocyte maintenance medium containing a vehicle control, a positive inducer (e.g., rifampicin for CYP3A4, Omeprazole for CYP1A2, Phenobarbital for CYP2B6, etc.), or test compounds.

Step 12: Return the plates to the incubator overnight.

Step 13: Change the medium with test compounds or a positive inducer once daily for three consecutive days.

Day 5

Step 1: Evaluate CYP enzymatic activity using preferred methods.

Step 2: Isolate total RNA if gene expression analysis is needed.

Day 1	Thaw and plate cells Corning Matrigel® matrix overlay (optional)
Day 2	Change to induction medium
Day 3	Change to induction medium
Day 4	Change to induction medium
Day 5	Evaluate CYP enzymatic activity Isolate RNA for gene expression analysis

Table 2. Recommended schedules for induction study

Procedure for hepatotoxicity studies

Day 2

Step 1: Remove the culture medium from each well and replace with pre-warmed Corning hepatocyte maintenance medium that contains test compounds and vehicle controls.

Step 2: Return the plates to the incubator overnight.

Day 3

Perform cytotoxicity assay using preferred methods.

Warranty/Disclaimer: Unless otherwise specified, all products are for research use only. Not intended for use in diagnostic or therapeutic procedures. Not for use in humans. Corning Life Sciences makes no claims regarding the performance of these products for clinical or diagnostic applications.

Mediatech, Inc.
A Corning Subsidiary
9345 Discovery Boulevard
Manassas, VA 20109
t 800.235.5476
t 703.471.5955
f 703.467.9851

[www.corning.com/
lifesciences/media](http://www.corning.com/lifesciences/media)

For additional product or technical information, please visit www.corning.com/lifesciences/media or call 1.800.235.5476.

CORNING | **FALCON**® | **AXYGEN** | **GOSELIN**™ | **PYREX**®

For a listing of trademarks, visit us at www.corning.com/lifesciences/trademarks. All other trademarks in this document are the property of their respective owners.